omak
www.omak.com.tr

63X3X

KULLANIM KILAVUZU
INSTRUCTION MANUAL

01.05.2016

63X3X KULLANIM KILAVUZU

**OMAK MAKİNA SANAYİİ ve TİCARET LİMİTED ŞİRKETİ
DR. MEDİHA ELDEM SOKAK 69/3 KOCATEPE /ANKARA
TEL : 312 433 31 33 (PBX) FAKS : 312 434 05 22
E-POSTA : omak@omak.com.tr WEB : www.omak.com.tr**

**omak**www.omak.com.tr

MENÜ AÇIKLAMALARI

SICAKLIK :

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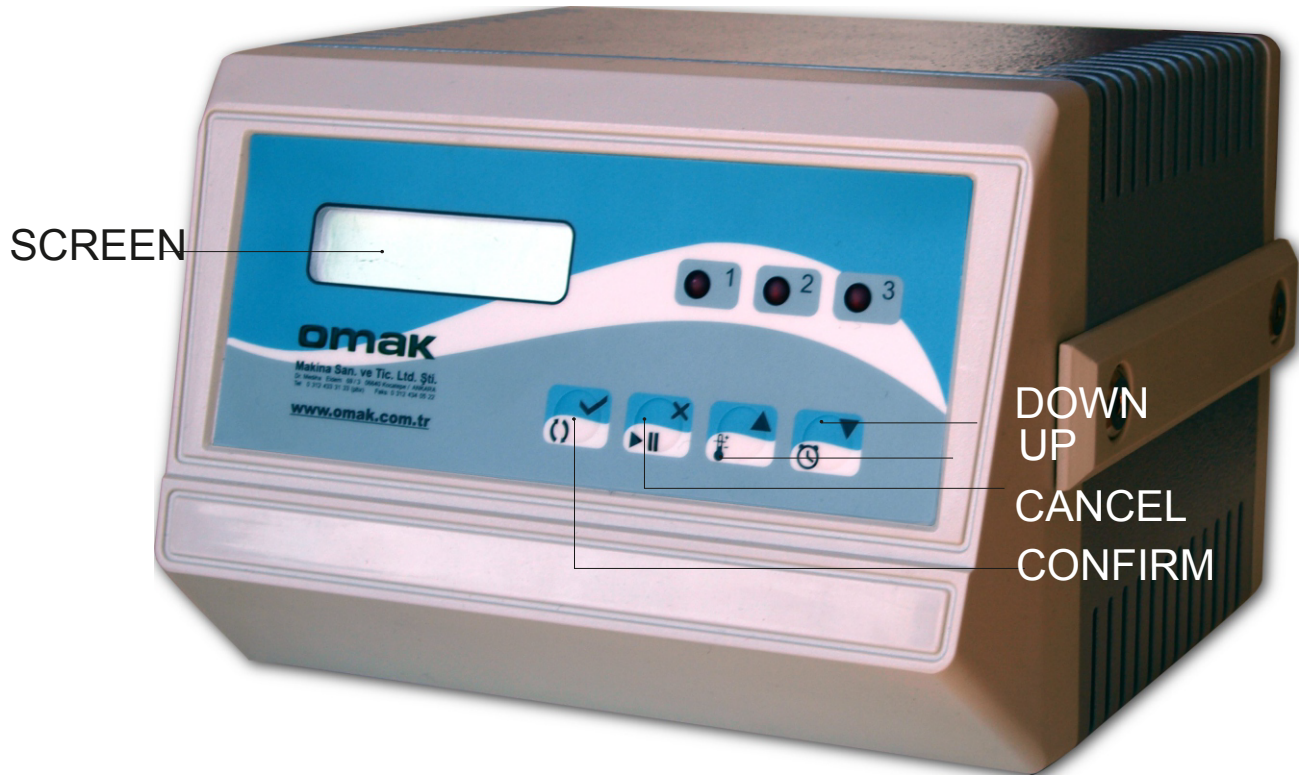
MENU INSTRUCTIONS

SICAKLIK (TEMPERATURE)

- S-SET_0** : Common temperature value. When the temperature value is defined from this section, the temperature values of the upper, middle and lower parts are automatically equalized to the value, set here.
- S-SET_U** : Used to define a value only for the upper section of the heater.
- S-SET_O** : Used to define a value only for the middle section of the heater.
- S-SET_A** : Used to define a separate value to the bottom section of the heater.
If all three sections are not wanted to be adjusted separately, they can be adjusted to the same temperature by using the first menu (S-SET_0),
- S-UST** : May assign a maximum limit temperature value for the heaters for safety.
For example, for a material that may be damaged at 80 C, S-UST value can be set to 65C therefore preventing any increase in temperature further.
- S-ALT** : May assign a minimum limit temperature value for the heaters for safety.
For example, for a material that may be melted above 40C, S-ALT value can be set to 40 C therefore preventing anybody to set a value below 40 C inadvertently.
- S-HSTZ** : In 63X3X; zero crossing semi conductor power control components are used as in the other devices of our range. Therefore, it does not produce any other noise. However, if the user needs a difference in temperature between the engagement and the disengagement, the user may define this hysteresis deficit to the device by S- HSTZ.

ÇIKIŞ (OUTPUT)

- C-UST** : the on / off setting of the upper section is made from this menu.
- C-ORTA** : the on / off setting of the middle section is made from this menu.
- C-ALT** : the on / off setting of the bottom section is made from this menu.

MENU EXPLANATIONS

ZAMAN (TIME)

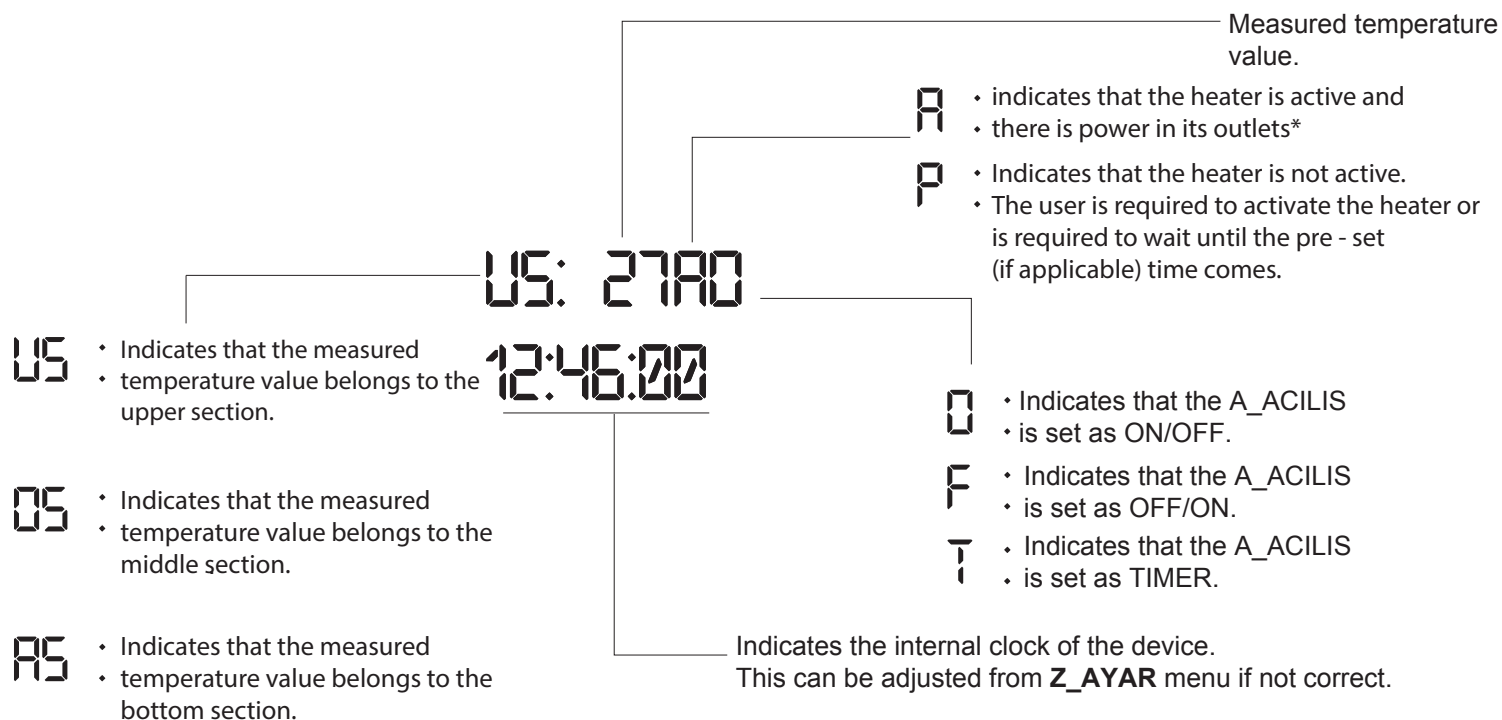
- Z-ACILIS** : The turn on time in automatic operation is set from this menu.
- Z_C.SURE** : We indicate how long a timer program, in which the turn on time is pre-set, will run from this menu.
- Z_AYAR** : Current time input is made to the device from this menu.

AYARLAR (SETTINGS)

- A_ACILIS** : 63XXX series present 3 different start up opportunities. The user may want to start the device by his own instead of auto start when the device is plugged in. In such a case, he should set the A_ACILIS setting to OFF/ON. If he wants the device heated up when the electricity is fed to the device, he should set it to ON/OFF. If he wants the device to wait until the preset starting time for heating up, he should set it as TIMER.
- A_KILIT** : The user may request the protection of the 63X3X device settings against any intervention. For this, you can lock the device from A_KILIT menu. The user cannot access to the menu in a locked device. In order to unlock, the energy should be fed to the device when the cancel key is pressed.
- A_VERS.** : This is the section, from where the version information is displayed.


SCREEN SHOTS

MAIN SCREEN SHOTS

It is at the user's discretion to halt the operation of the device or resuming the operation of a device in standby mode when the device is on/off or off / on mode. This option can be used by pressing cancel key on the main menu.

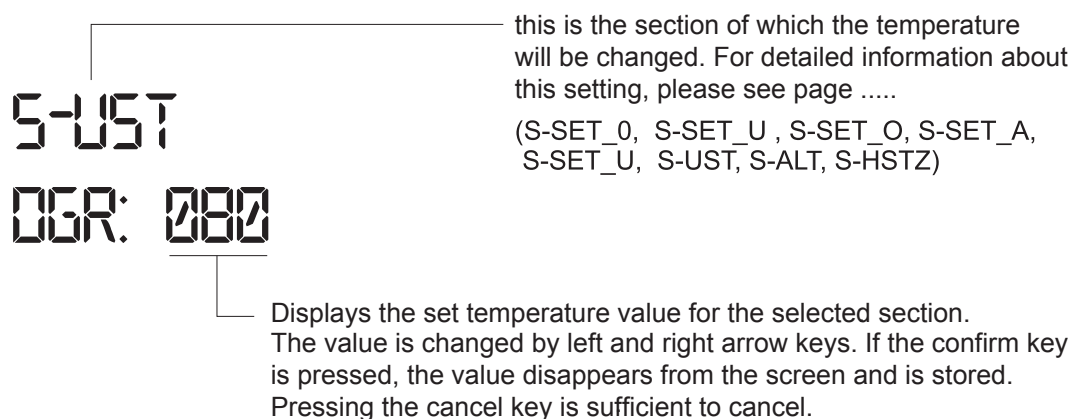TEMPERATURE SETTINGS MENU

OUTLET SETTINGS MENU

C-ALT
ACK

Shows the section from which the outlet adjustment shall be made. Those sections may be one of the following. For detailed information about this setting, please see page
(C-UST, C-ORTA, C-ALT)

You can power on and power off the selected section of the device by up and down arrow keys and you can store this setting with Accept key. After storing the settings, ON or OFF will be disappeared from the screen. The Cancel Key, as in other menus, exits the current menu without completing the storing the settings.

TIME SETTINGS MENU

Z-ACILIS
00:XX

Z-ACILIS, Z-C.SURE and Z-AYAR menus are stored by the same method. The clock is adjusted when XX characters are on the right side. After storing the hour value with accept key, digits appear in place of XX characters. After the hour characters appear, the minute value should be set with up / down arrow keys.

On the right side, XX characters appear on the right side when the hour value is set.

Z-ACILIS
00:00

After the hour value, set by the accept key, minute appears on the right side instead of XX characters.

The minute value is changed by up and down arrow keys and the value is stored by the accept key.

SETTINGS MENU

A-ACILIS
ON/OFF

Start settings section determines the operation mode in the device start up. You can set it as **ON / OFF**, **OFF / ON** or **TIMER** by the up / down arrow keys.

The selected value will appear in the bottom line.

A-KILIT
ACIK

It is possible to lock the device in order to prevent changing the device settings. After the device lock, a user cannot access to the menus. In order to unlock, first the cancel key is pressed and power is fed to the device. In the screen, **KILIT ACIK** will appear.

The name of the process to be undertaken is written in the bottom line.

OMAK MAKİNA SANAYİİ ve TİCARET LİMİTED ŞİRKETİ
DR. MEDİHA ELDEM SOKAK 69/3 KOCATEPE /ANKARA
TEL : 312 433 31 33 (PBX) FAKS : 312 434 05 22
E-POSTA : omak@omak.com.tr WEB : www.omak.com.tr

omak

www.omak.com.tr